**GPS SERIES USER MANUAL/  
BENUTZERHANDBUCH  
GPS-REIHE/MANUEL DE  
L'UTILISATEUR SÉRIE GPS/  
MANUAL DEL USUARIO  
DE LA SERIE GPS/ MANUALE  
UTENTE SERIE GPS**

**Laboratory Refrigerators and Freezers**

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**-20°, + 4°**

## Manufacturer's Responsibility

The Manufacturer is responsible as to safety and correct use of these refrigerators only if:

- the operations of setting, change, and repair must be performed by authorised personnel.
- the electrical installation has been carried out in conformity with the instructions.
- the device is used in accordance with this manual.

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### **SPECIAL WARNINGS:**

Not to be used in the presence of explosive gases or mixtures and / or close to sources with high magnetic or electric fields. Not intended for flammable material storage.

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## **COPYRIGHT & NOTES**

All rights reserved. No part of this document is to be reproduced, stored, or transmitted in any form without the prior written permission of the manufacturer. Despite every care being adopted during preparation of the document. The manufacturer shall not assume any responsibility for atrocities or negligence. Nor shall it assume any responsibility for damages caused by use of the information contained herein. These same rights are reserved for any person(s) or company involved in the creation and production of this manual. All trademarks mentioned in this manual are property of the respective companies.

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## **Index**

	<b><u>Products applicable to this manual</u></b>	<b>6</b>
<b>1</b>	<b>Standards and general warnings</b>	<b>6</b>
1.1	Testing and intended use	
1.2	Introduction	
1.3	Product description	
1.4	General safety regulations	
1.5	Client's responsibilities	
1.6	Client service requests	
1.7	Ordering spare parts	
1.8	Product configuration	
1.9	Materials and refrigerants	
<b>2</b>	<b>Installation</b>	<b>24</b>
2.1	Transport and handling	
2.2	Positioning	
2.3	Wiring and electrical hook-up	
2.4	Set-up operations	
2.5	Re-installation	
2.6	Scrapping and disposal	
<b>3</b>	<b>Safety</b>	<b>35</b>
3.1	Safety and accident prevention	
3.2	Safety dataplates and guards	
3.3	Operating limits	
3.4	Conditions of storage	
<b>4</b>	<b>User Instructions</b>	
<b>4</b>	<b>Operation</b>	<b>40</b>
4.1	General warnings	
4.2	Controller general description	
4.3	Regulation	
4.4	Defrost	
4.5	Control of evaporator fans	
4.6	Front panel commands	

- 4.6.1 Use of LEDs
- 4.7 Alarm signaling
  - 4.7.1 Alarm recovery
- 4.8 Technical data
- 4.9 Connections
- 4.10 Default setting values

**5 Routine maintenance 76**

---

- 5.1 Cleaning the refrigerator
  - 5.1.1 Cleaning the interior and exterior of the equipment
  - 5.1.2 Cleaning the condenser
- 5.2 Precautionary measures for prolonged disuse

**6 Troubleshooting 83**

---

- 6.1 Troubleshooting

**7 Warranty 89**

---

# 1 STANDARDS AND GENERAL WARNINGS / NORMEN UND ALLGEMEINE WARNHINWEISE / NORMES ET AVERTISSEMENTS GÉNÉRAUX / NORMAS Y ADVERTENCIAS GENERALES / AVVISI STANDARD E GENERALI

## PRODUCTS APPLICABLE TO THIS MANUAL / AUF DIESES HANDBUCH ANWENDBARE PRODUKTE / PRODUITS APPLICABLES À CE MANUEL / PRODUCTOS APLICABLES A ESTE MANUAL / PRODOTTI TRATTATI IN QUESTO MANUALE

The present manual is exclusively valid and applicable to the following products series / Die Geltung und Anwendbarkeit dieses Handbuchs erstreckt sich ausschließlich auf folgende Produktreihen / Le présent manuel est exclusivement valide et applicable aux séries de produits suivantes / Este manual solo es válido y aplicable a las siguientes series de productos / Il presente manuale vale e si riferisce esclusivamente alle seguenti serie di prodotti:

### GPS Series Laboratory refrigerators / Laborkühlschränke der GPS-Reihe / Réfrigérateurs de laboratoire série GPS / Frigoríficos de laboratorio de la serie GPS / Frigoriferi da laboratorio serie GPS

**Adjustable temperature control range:** lowest T = 2 °C, highest T = +16°C

**Operating temperature:** +4°C to 6°C

**Factory pre-set to:** +4°C

**Models:** R400-SAEV/W, R400-GAEV/W, R700-SAEV/W, R700-GAEV/W, R14X-SAEV/W, R14X-GAEV/W

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**Modelle:** R400-SAEV/W, R400-GAEV/W, R700-SAEV/W, R700-GAEV/W, R14X-SAEV/W, R14X-GAEV/W

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**Preimpostata in fabbrica a:** +4°C

**Modelli:** R400-SAEV/W, R400-GAEV/W, R700-SAEV/W, R700-GAEV/W, R14X-SAEV/W, R14X-GAEV/W

# 1 STANDARDS AND GENERAL WARNINGS / NORMEN UND ALLGEMEINE WARNHINWEISE / NORMES ET AVERTISSEMENTS GÉNÉRAUX / NORMAS Y ADVERTENCIAS GENERALES / AVVISI STANDARD E GENERALI

**GPS Series Laboratory freezers / Laborgefrierschränke der GPS-Reihe / Congélateurs de laboratoire série GPS / Congeladores de laboratorio de la serie GPS / Congelatori da laboratorio serie GPS**

**Adjustable temperature control range:** lowest T = -25 °C, highest T = -10°C

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All relevant data referring to these products can be found on the data label visible on the rear part of the cabinet. Here is an example of the label:

<b>Mod:</b> R400-GAEV	
<b>Serial Number:</b> 11POP 004349	
<b>Month, Year:</b> Jul/2011	
<b>Voltage:</b> 220-230V-1~	
<b>Frequency:</b> 50 Hz	
<b>Power:</b> 430 W	
<b>Refrigerant Gas Type:</b> R404A	
<b>Gas Q.ty:</b> 0.300 kg	
<b>Brut Vol:</b> 1050 l	
<b>Net Vol:</b> 700 l	
<b>Group No:</b> 1	
	
R400-GAEV 11POP 004349	
<b>MADE IN ITALY</b>	

# 1 STANDARDS AND GENERAL WARNINGS / NORMEN UND ALLGEMEINE WARNHINWEISE / NORMES ET AVERTISSEMENTS GÉNÉRAUX / NORMAS Y ADVERTENCIAS GENERALES / AVVISI STANDARD E GENERALI

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This equipment is tested in compliance with established regulations and then shipped ready for use. This equipment is intended for general purpose laboratory cold storage.

*"If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired."*

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## 1.2 INTRODUCTION / EINFÜHRUNG / INTRODUCTION / PRESENTACIÓN / INTRODUZIONE

This manual provides all instructions required for the correct use of the equipment and to keep it in optimal condition. It also contains important user safety information.

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The following professional roles are explained in order to define individual responsibilities:

**Installer:** a qualified technician who installs the equipment in accordance with these instructions.

**User:** the person who, after having read this manual carefully, uses the equipment in accordance with the intended specification of use described in this manual.

User's responsibilities:

- ensure that the product is kept at suitable temperatures in an ambient environment less than +32° C;
- be aware of the regulations governing the conservation of products to refrigerate and to observe any whatsoever hygiene indications that may be applicable.

# 1 STANDARDS AND GENERAL WARNINGS / NORMEN UND ALLGEMEINE WARNHINWEISE / NORMES ET AVERTISSEMENTS GÉNÉRAUX / NORMAS Y ADVERTENCIAS GENERALES / AVVISI STANDARD E GENERALI

The user is obliged to carefully read the manual and refer to its information at all times. Particular attention must be paid to safety warnings (refer to Section 1.4).

**Routine maintenance technician:** qualified operator able to perform routine maintenance of the equipment by following the instructions in this manual (can be the customer too).

**Service engineer:** qualified technician, authorised by the manufacturer to perform extraordinary maintenance of the equipment .

The symbol  appears at certain points in the manual to draw the reader's attention to important safety information.

The manufacturer declines any responsibility in case of improper use of the equipment deviating from the reasonably construed intended use, and for all operations carried out that are not in compliance with the instructions reported in the manual.

This manual must be stored in an accessible and known place for all operators (installer, user, routine maintenance technician, service engineer).

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The equipment comprises a single body with panelling in various materials and insulation in expanded polyurethane foam density. The equipment instruments are located on the front panel where the electrical wiring is housed. The motor unit and the condenser unit are housed on the top of body. The interior parts are fitted with suitable supports for shelves. The doors are fitted with an automatic return device and magnetic seal elements. During the design and construction stage all measures have been adopted to implement total safety including radius interior corners, funnel-shaped base panel to convey condensate to exterior, no rough surfaces, fixed guards protecting moving or potentially dangerous parts.

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Read this manual carefully and follow the prescriptions contained herein.

The user assumes full responsibility in case of operations carried out without observing the instructions in the manual.

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**Do not use this product with flammable gases or solvents that burn easily.**

Primary general safety regulations:

- Do not touch the unit with wet hands and/or feet
- Do not use the equipment with bare feet
- Do not insert screwdrivers or other pointed objects between guards or moving parts of the equipment
- Do not pull the power cord to disconnect the equipment from the electrical mains
- Make sure that the equipment is not used by unsuitably qualified persons
- Before performing any cleaning or maintenance on the equipment disconnect it from the electrical mains by switching off the main switch and extracting the plug
- In case of faults or malfunctions, switch off the equipment and do not attempt to repair it by yourself as doing so may void the warranty. All service and repair operations must be performed exclusively by a suitable authorized engineer.

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The customer is required to:

- Execute the electrical connection of the equipment
- Prepare the place of installation
- Provide consumable materials for cleaning
- Perform routine maintenance

In the case of power failures or malfunctions do not open the doors, in order to maintain as much as possible the internal temperature. If the problem persists for more than a few hours, move the contents to a more suitable place.

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For all technical problems and any requests for technical service, refer exclusively to the contact information on the last page of this manual or your local dealer.

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Orders of spare parts should be made by consulting the part reference code and the serial number of your unit. Consult your dealer.

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The unit is designed solely for the preservation of laboratory products, which require various controls and warning in case of sudden alteration of temperature.

# 1 STANDARDS AND GENERAL WARNINGS / NORMEN UND ALLGEMEINE WARNHINWEISE / NORMES ET AVERTISSEMENTS GÉNÉRAUX / NORMAS Y ADVERTENCIAS GENERALES / AVVISI STANDARD E GENERALI

**! PRODUCTS MUST BE STORED IN ORDER TO ENSURE EFFICIENT AIR CIRCULATION INSIDE THE UNIT AND SHALL NOT COME OUT OF THE SHELF PERIMETER.**

All uses outside of manufacturer's intended use in section 1.1 shall be construed as "improper use" for which the manufacturer declines all responsibility.

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Materials in contact or potentially in contact with products are in compliance with the relevant directives. The equipment is designed and built so that contact parts can be cleaned before each use. The refrigerants utilized comply with established regulations.

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**!** The equipment must be transported and handled exclusively in upright position, in observance of the instructions printed on the packing.

This precaution is necessary to avoid contamination of the refrigerant circuit with compressor lube oil with resulting valve and heat exchanger coil failure and problems starting the electric motor.

The manufacturer is not responsible for any problems due to transport executed in conditions other than those specified herewith.

The equipment is secured to a Styrofoam base and wrapped in a plastic film.

The equipment must be handled using a fork lift truck or a pallet truck with suitable forks (fork length at least equal to 2/3 length of unit).

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Incorrect positioning can cause damage to the equipment and generate hazardous conditions for personnel. The installer must therefore observe the following general regulations:

- Make sure you maintain a minimum of 10 cm. clearance from the walls and the ceiling
- The room must be well ventilated
- Keep well away from sources of heat
- Avoid direct sunlight
- Remove packing material
- Remove accessories from inside the unit
- Styrofoam or Wood base removal: using a hammer, tilt the cabinet to one side and loosen the two thread-forming screws, drag the cabinet from the back side holding the base still until the four feet have gone out from the containing holes, slightly tilt the cabinet backward and take the base away pulling it from the front side.
-  Use gloves when handling Styrofoam or wooden packing materials and the wooden base to protect the hands from splinters.
- Position the equipment with the help of a level. Remove the protective PVC film from the external surfaces of the unit
- Position the shelf runners in the holes in the uprights
- Insert the shelves in the runners
- Insert the condensate collection tray, if needed, in the runners located beneath the unit

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The electrical plant and electrical hook-up operations must be performed by a qualified electrician. For safety reasons adhere to the following indications:

- Check that the electrical plant is suitably sized for the absorbed power of the unit
- If the electrical socket and the plug on the equipment power cord are incompatible, call technical service (see last page) or your local distributor.
- Do not use reductions or multi-way adapters (Fig.1)
-  It is important to connect the equipment correctly to an efficient earth system executed in compliance with the relevant legislation.
- The equipment must be positioned so that plug can be easily reached (Fig. 1)

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To avoid errors and accidents, perform a series of checks for possible damage sustained during transport, installation and hook-up operations before starting up the unit.

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#### *PRELIMINARY CHECKS*

- Check the condition of the power cord (no cut or chaffing)
- Check that the door hinges and shelf support are stable
- Check the door seals and shelves are not damaged (broken or scratched) and that the door closes and seals properly
- Make sure pipelines, unions are in perfect condition

#### *INDICATIONS FOR OPTIMAL DUTY*

- Do not block the motor compartment air vents
- Do not lay objects on the top of the equipment
- Before storing products wait until they are cold
- Arrange the products on suitable shelves or in containers. Do not place products directly on the base or against the walls, doors or fixed guards of the unit
- Make sure doors are kept closed
- Keep the defrost water drain outlet clear
- Limit the frequency and duration of opening; each time the door is opened the internal temperature will alter
- Load products at ambient temperature gradually to allow correct refrigeration
- Perform routine maintenance regularly (see sect. 5)

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Observe the following procedure:

- Switch off the equipment from the main switch
- Disconnect the power cord from the electrical outlet
- Handle the equipment in accordance with the instructions in Section 2.1
- Follow the instructions in Section 2.2 for positioning and hook-ups in the new location

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These units may contain materials, which at the end of the working life of the apparatus, must be disposed at one of the recycling centres nominated by your Local National Health Department or as specified by the law in force. Scrapping and disposal of the equipment must be carried out in full observance of established legislation in your country.

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In particular, the apparatus may contain the following materials:

- Iron
- Copper
- Aluminium
- Non-biodegradable plastics
- Fibre glass for printed circuits
- Ferrite
- Batteries
- CFC-free refrigeration gas
- Electrical and electronic equipment (WEEE)

The manufacturer shall not be chargeable for any disposal of the apparatus at the end of its working life.

In line with EU Directive 2002/96/EC for waste electrical and electronic equipment (WEEE), this electrical product must not be disposed of as unsorted municipal waste. Please dispose of this product by returning it to your local municipal collection point for recycling.

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Refrigerators and Freezers are designed to include several features to ensure the safety and the protection of the operators and for the material stored in the cabinets.

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The following list describes the protections adopted against mechanical risks:

- **stability:** the equipment is designed and built so that even with the shelves fully extended in the intended conditions of operation, the unit will remain stable, so that it can be used without risk of tipping, falling or sudden movement.
- **surface, edges, corners:** accessible parts of the equipment have no sharp corners, sharp edges or rough surface that could cause injuries.
- **moving parts:** moving parts of the unit are designed, built and configured to avoid risks. Moving parts are protected by fixed guards to prevent accidental contact that could result in injury.

Measures adopted for protection against additional risks:

- **electrical power:** the unit is designed, built and fitted out with the aim of preventing the risk of electric shock in compliance with established safety legislation
- **noise:** the equipment is designed and built to reduce risks related to the emission of airborne noise to the minimum level.

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**It is strictly forbidden:**

- To tamper or remove the evaporator cover that protects the user from the risk of being cut on the heat exchanger fins;
- To remove the data plate fixed to the inside edge of the motor housing showing technical specification and earth connection warning;
- To remove the data plates on the evaporator unit cover near the electrical wiring inside the motor housing which warn the user to disconnect electrical power before working on unit;
- To remove the data tag fixed to the power cord showing the type of power supply.

The manufacturer declines all responsibility for safety of the equipment if the above recommendations are not observed.

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This equipment is designed and built to work with ambient temperatures between +10°C and +32°C. Units are designed and tested for temperate countries.

If the ambient conditions are different it will not be possible to achieve the top performance levels specified by the manufacturer.

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The refrigerator if not used must be stored indoors in a temperature not exceeding +40 ° C or below 0 ° C.

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### **4 OPERATION / BETRIEB / FONCTIONNEMENT / FUNCIONAMIENTO / FUNZIONAMENTO**

Before switching on the unit, check that the electrical connections have been made correctly and above all, that the ground connection is available and working properly. Perform preliminary cleaning of the unit as described at Section 5.

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##### **Please read before using this manual**

- This manual is part of the product and should be kept near the instrument for easy and quick reference.
- Digital controller with defrost and fans management shall not be used for purposes different from those described hereunder. It cannot be used as a safety device.
- Check the application limits before proceeding.

##### **Safety precautions**

- Check the supply voltage is correct before connecting the instrument.
- Do not expose to water or moisture: use the controller only within the operating limits avoiding sudden temperature changes with high atmospheric humidity to prevent formation of condensation
- Warning: disconnect all electrical connections before any kind of maintenance.
- In case of failure or faulty operation contact technical service or your local distributor.
- Consider the maximum current which can be applied to each relay
- Ensure that the wires for probes, loads and the power supply are separated and far enough from each other, without crossing or intertwining.

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The controller, format 32 x 74 x 60 mm, is microprocessor based controller, suitable for applications on medium or low temperature ventilated refrigerating units. It has three relay outputs to control compressor, fan, and defrost which can be either electrical or reverse cycle (hot gas). The device is also provided with 2 NTC or PTC probe inputs, the first one for temperature control and the second one to be located onto the evaporator, to control the defrost termination temperature and to managed the fan, the third one, optional, to connect to the HOT KEY terminals to signal the condenser temperature alarm or to display a temperature. The digital input can operate as fourth temperature probe.

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
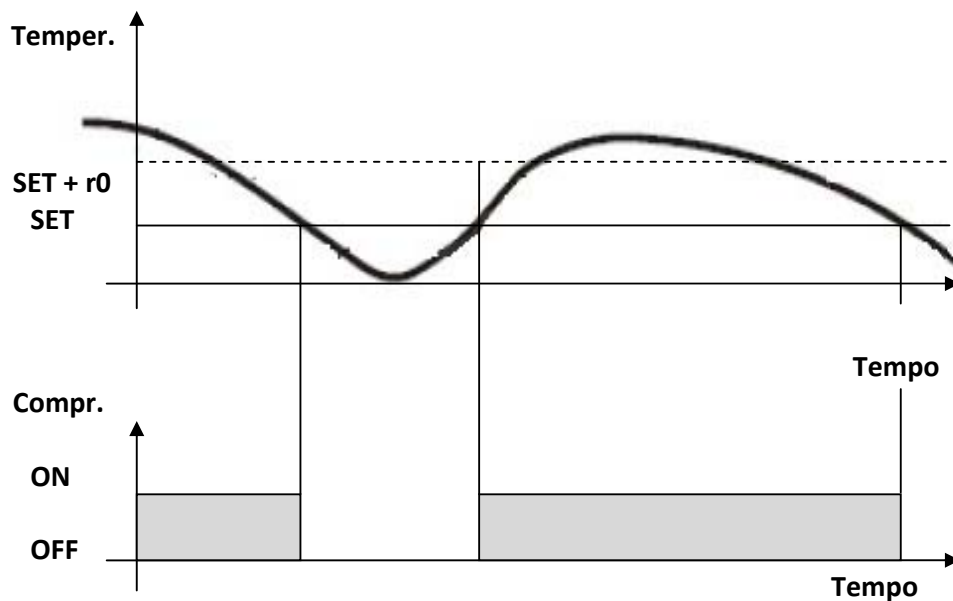
The regulation is performed according to the temperature measured by the thermostat probe with a positive differential from the set point: if the temperature increases and reaches set point plus differential the compressor is started and then turned off when the temperature reaches the set point value again.

In case of fault in the thermostat probe the start and stop of the compressor are timed through parameters "Cy" and "Cn".

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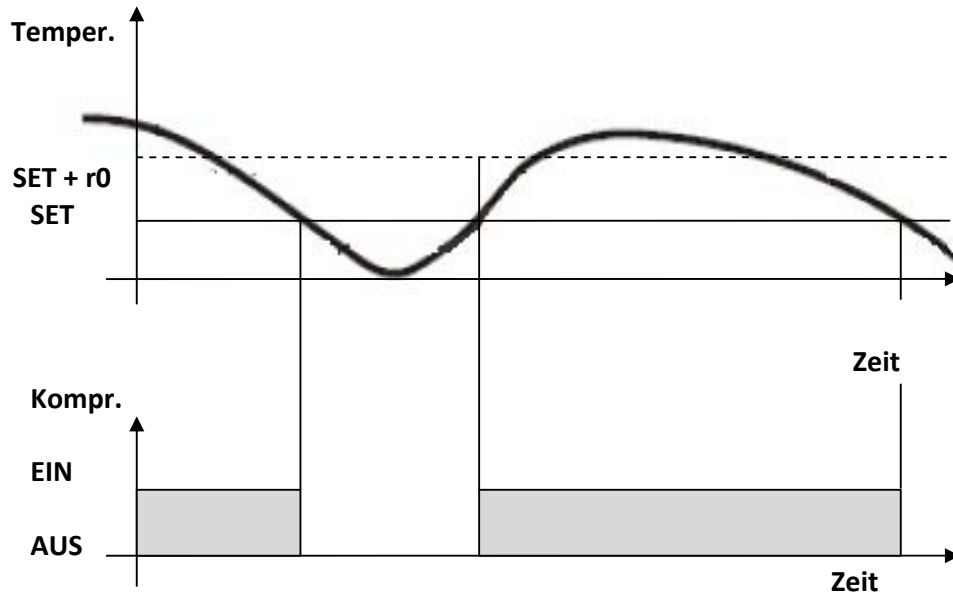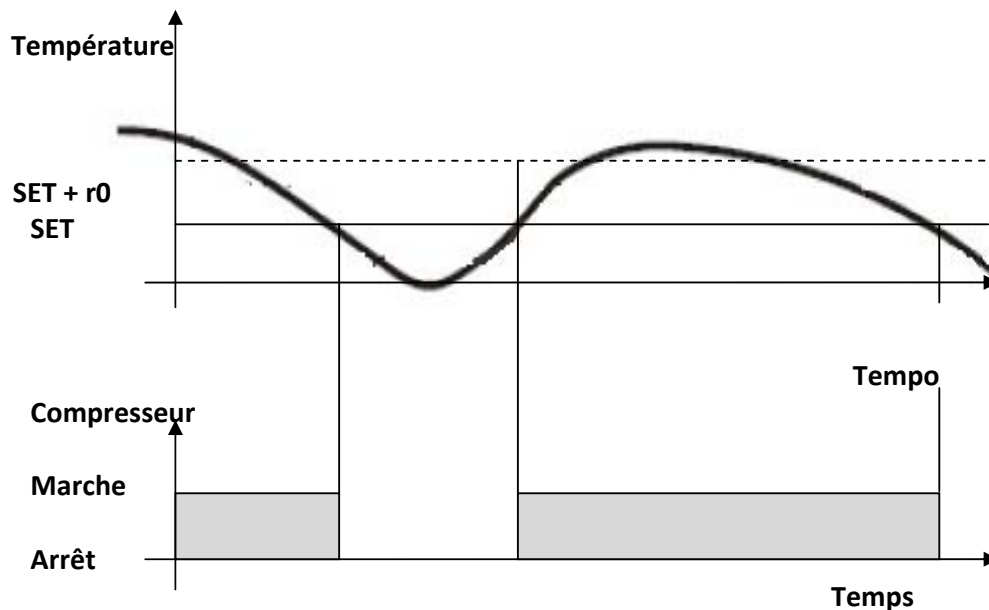
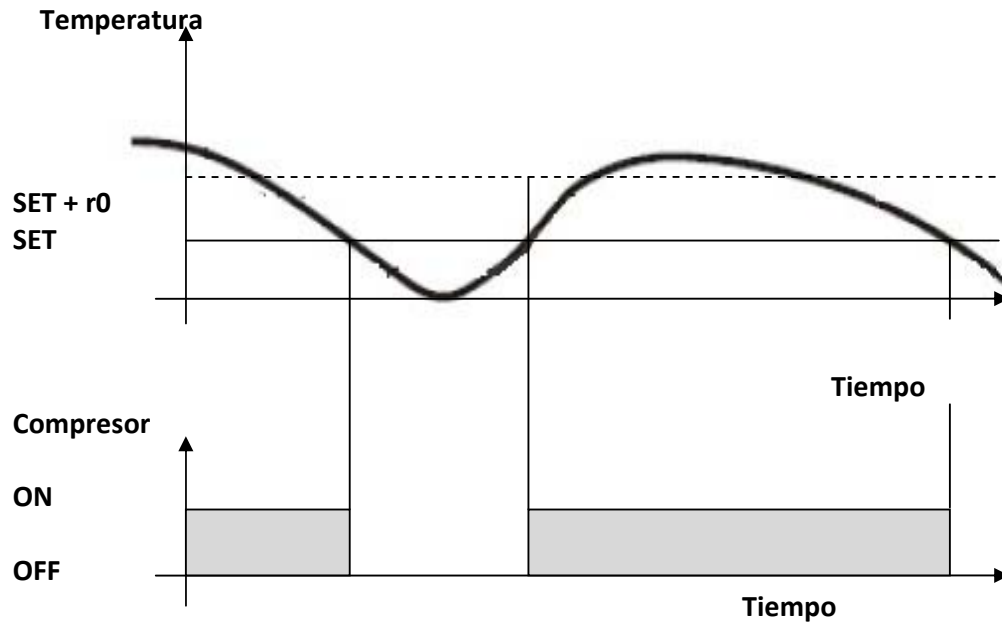
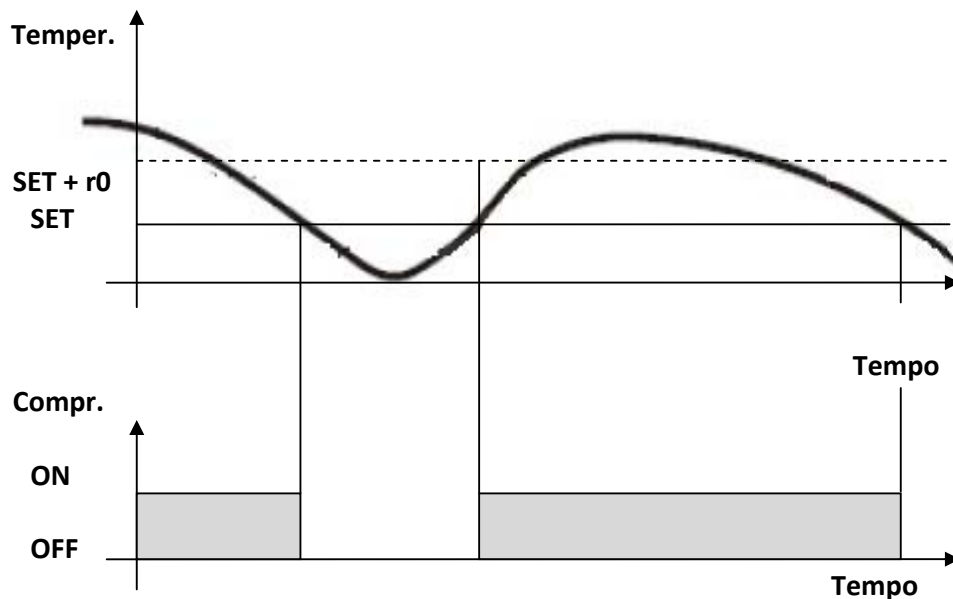
#### 4.4 DEFROST / ABTAUEN / DÉGIVRAGE / DESCONGELACIÓN / SBRINAMENTO

The factory default setting of the defrost parameter is tdf=in (hot gas defrost) for both refrigerators and freezers

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The fan control mode is selected by means of the “FnC” parameter: The factory default setting is o\_n.

- **FnC = o\_n** fans will run even if the compressor is off, and not run during defrost; After defrost, there is a timed fan delay allowing for drip time, set by means of the “Fnd” parameter.

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**SET**: To display target set point; in programming mode it selects a parameter or confirm an operation.

❄ : **(DEF)** To start a manual defrost

▲ **(UP)**: To see the max. Stored temperature; in programming mode it browses the parameter codes or increases the displayed value.

▼ **(DOWN)**: To see the min. Stored temperature; in programming mode it browses the parameter codes or decreases the displayed value.

⏻  
💡 To switch the instrument off, if onF=OFF.  
Not enabled.

##### KEY COMBINATIONS:

▲ + ▼ To lock & unlock the keyboard.

SET + ▼ To enter in programming mode.

SET + ▲ To return to the room temperature display.

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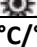
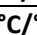
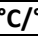
Each LED function is described in the following table.

LED	MODE	FUNCTION
	ON	Compressor enabled
	FLASHING	Anti-short cycle delay enabled
	ON	Defrost enabled
	FLASHING	Drip time in progress
	ON	Fans enabled
	FLASHING	Fans delay after defrost in progress
	ON	An alarm is occurring
	ON	Continuous cycle is running
	ON	Energy saving enabled
°C/°F	ON	Measurement unit
°C/°F	FLASHING	Programming phase

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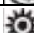
### How to see the min temperature

1. Press and release the down key.
2. The "Lo" message will be displayed followed by the minimum temperature recorded.
3. By pressing the down key again or by waiting 5s the normal display will be restored.

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### How to see the max temperature

1. Press and release the up key.
2. The "Hi" message will be displayed followed by the maximum temperature recorded.
3. By pressing the up key again or by waiting 5s the normal display will be restored.

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### **How to reset the max and min temperature recorded**

1. Hold press the SET key for more than 3s, while the max. or min temperature is displayed. (rSt message will be displayed)
2. To confirm the operation the "rSt" message starts blinking and the normal temperature will be displayed.

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### **How to see the set point**

1. Push and immediately release the **SET** key, the set point will be showed;
2. Push and immediately release the **SET** key or wait about 5s to return to normal visualization.

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#### **How to change the set point**

**Units are factory pre-set to the correct temperature. Changing the set point is not recommended.**

1. Push the SET key for more than 2 seconds to change the Set point value;
2. The value of the set point will be displayed and the "°C" or "°F" LED starts blinking;
3. To change the value, press the ▲ or ▼ arrows.
4. To store the new set point value in memory, push the SET key again or wait 10 seconds.

#### **How to start a manual defrost**

Push the **DEF** key for more than 2 seconds and a manual defrost will start

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#### How to change a parameter value

To change the parameter's value operate as follows:

1. Enter the Programming mode by pressing the **SET+ ▼** keys for 3s (“°C” or “°F” LED starts blinking).
2. Select the required parameter. Press the “**SET**” key to display its value
3. Use ▲ or ▼ to change its value.
4. Press “**SET**” to store the new value and move to the following parameter.

**To exit:** Press **SET+ ▲** or wait 15s without pressing a key.

**NOTE:** the set value is stored even when the procedure is exited by waiting the time-out to expire.

**Parameter changes should only be performed by qualified Service personnel. Refer to the Service section at the back of this manual for controller details.**

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Message	Cause	Outputs
"P1"	Room probe failure	Compressor output acc. To par. "Con" and "COF"
"P2"	Evaporator probe failure	Defrost end is timed
"P3"	Third probe failure	Outputs unchanged
"P4"	Fourth probe failure	Outputs unchanged
"HA"	Maximum temperature alarm	Outputs unchanged
"LA"	Minimum temperature alarm	Outputs unchanged
"HA2"	Condenser high temperature	It depends on the "Ac2" parameter
"LA2"	Condenser low temperature	It depends on the "bLL" parameter
"dA"	Door open	Compressor and fans restarts
"EA"	External alarm	Output unchanged
"CA"	Serious external alarm(i1F=bAL)	All outputs OFF
"CA"	Pressure switch alarm(i1F=PAL)	All outputs OFF

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Probe alarms "P1", "P2", "P3" and "P4" start some seconds after the fault in the related probe; they automatically stop some seconds after the probe restarts normal operation. Check connections before replacing the probe.

Temperature alarms "HA", "LA", "HA2" and "LA2" automatically stop as soon as the temperature returns to normal values.

Alarms "EA" and "CA" (with i1F=bAL) recover as soon as the digital input is disabled.

Alarm "CA" (with i1F=PAL) recovers only by **switching off and on** the instrument.

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### **Basic safety regulations**

We summarise the elementary safety regulations to ensure that the user or maintenance technician can perform the work in conditions of total safety:

- do not touch the unit with wet hands and/or feet;
- do not use the equipment with bare feet;
- do not insert screwdrivers or other pointed objects between guards or moving parts of the equipment;
- do not pull the power cord to disconnect the equipment from the electrical mains;
- before performing any cleaning or maintenance on the equipment disconnect it from the electrical mains by switching off the main switch and extracting the plug.

### **Prohibited: removal of guards and safety devices**

It is strictly forbidden to remove guards or safety devices when performing routine maintenance work. The manufacturer disclaims all liability that may arise if this regulation is not observed.

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### **Emergency measures in case of fire**

- Disconnect the unit from the electrical power socket.
- Do not use water to extinguish the fire.
- Use powder or foam extinguishers.

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The unit is designed to preserve laboratory products so it is important to keep it clean. The equipment is thoroughly cleaned in our factory before delivery. We recommend, however, that you clean the interior of the equipment before use. Before cleaning the equipment, make sure the power cord is disconnected.

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The condenser is a heat exchanger. If it is dirty or clogged the air cannot pass freely into the same, it cannot do its job properly and the efficiency of the refrigeration system falls down proportionally. **FOR THOSE REASONS IT IS IMPORTANT TO KEEP CLEAN THE CONDENSER, TYPICALLY MONTHLY.**

#### **WARNING!!**

**Always switch off the unit and disconnect power cord before cleaning, it is dangerous to do it with power: fan may start suddenly at any time!**

Use a convenient ladder to reach the condenser. Use an air jet or an aspirator with a soft dry brush if necessary (Fig. 12) and remove any dust or fluff from the heat exchanger fins.

After cleaning, start the equipment.

**/!** During this operation use the following personal safety measures:

- safety glasses,
- respirator mask,
- gloves

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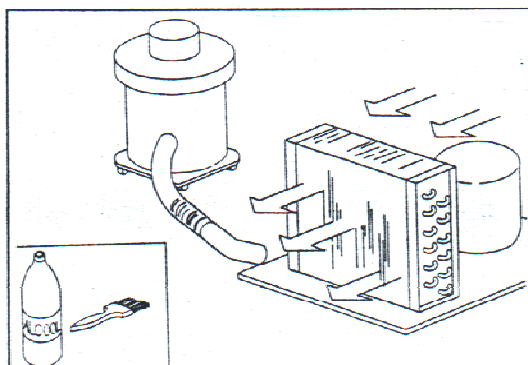
If the refrigerator should remain unused for more than 15 days please proceed as follows:

- Switch off the equipment and disconnect the power cord,
- Clean the interior of the cabinet, shelves, trays, runners and supports, paying special attention to critical areas such as articulations and magnetic sealing strips,
- Leave doors slightly open to prevent accumulation of residual humidity.

All maintenance work not described in the previous sections must be considered "Special Maintenance".

Special maintenance interventions and repairs are to be performed exclusively by specialised technicians authorised by the manufacturer.

The manufacturer declines all liability in the case of work performed by the user or unauthorised persons, or if non-original spare parts are fitted to the equipment.

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Below is a list of problems, with possible causes and recommended actions.

FAULT	POSSIBLE CAUSE	ACTION
The equipment does not switch on	Power switch is OFF	Move main switch ON
	Power failure	Check the plug, socket, fuses, electrical line
	Other	Contact service department
Refrigeration unit does not start	Set temperature has already been reached	Set new temperature
	Defrosting cycle in progress	Wait for the end of defrost cycle, switch off and then on again after few minutes
	Control panel fault	Contact service department
	Other	Contact service department
Refrigeration unit runs constantly but does not reach set temperature	Room temperature too high	Provide for better ventilation of the room and clean condenser
	Condenser dirty	Clean condenser
	Refrigerant needs recharging	Contact service department
	Fan of condenser does not work	Contact service department
	Door gasket is damaged	Check seals /make sure refrigerator contents are not obstructing the door
	Evaporator is coated with ice	Defrost manually
	Defrost valve is open	Contact service department
Water or ice deposits in the drip tray	Drain outlet is clogged	Clean drain and drain outlet
	Unit is not levelled	See heading 2.2

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## 7 WARRANTY / GARANTIE / GARANTIE / GARANTÍA / GARANZIA

### Thermo Scientific GPS Series Refrigerators and Freezers European Warranty Policy

---

---

#### **Thermo Scientific GPS Series Refrigerators • European Warranty • 12 Months Full Warranty Parts and Labor**

During the first twelve (12) months from shipment, Thermo Fisher Scientific Inc, through its authorized Dealer or service organizations, will at its option and expense repair or replace any part found to be non-conforming in material or workmanship. Thermo Fisher Scientific Inc reserves the right to use replacement parts, which are used or reconditioned. Replacement or repaired parts will be warranted for only the unexpired portion of the original warranty.

This warranty does not apply to damage caused by (i) accident, misuse, fire, flood or acts of God; (ii) failure to properly install, operate or maintain the products in accordance with the printed instructions provided, (iii) causes external to the products such as, but not limited to, power failure or electrical power surges, (iv) improper storage and handling of the products, (v) use of the products in combination with equipment or software not supplied by Thermo Fisher; or (vi) installation, maintenance, repair, service, relocation or alteration of the products by any person other than Thermo Fisher or its authorized representative. To obtain proper warranty service, you must contact the nearest authorized service center or Dealer. Thermo Fisher Scientific, Inc's own shipping records showing date of shipment shall be conclusive in establishing the warranty period. At Thermo Fisher's option, all non-conforming parts must be returned to Thermo Fisher postage paid and replacement parts are shipped FOB Thermo Fisher's location.

#### **Limitation of Liability**

THIS WARRANTY IS EXCLUSIVE AND IN LIEU OF ALL OTHER WARRANTIES, WHETHER WRITTEN, ORAL, OR IMPLIED. NO WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE SHALL APPLY. THERMO FISHER DOES NOT WARRANT THAT THE PRODUCTS ARE ERROR-FREE OR WILL ACCOMPLISH ANY PARTICULAR RESULT. THERMO FISHER SHALL NOT BE LIABLE FOR ANY INDIRECT OR CONSEQUENTIAL DAMAGES INCLUDING, WITHOUT LIMITATION, DAMAGES TO LOST PROFITS OR LOSS OF PRODUCTS.

**For technical or sales support please contact your local distributor or Thermo Scientific at the following numbers:**

**Europe:** Austria +43 1 801 40 0, Belgium +32 2 482 30 30, France +33 2 2803 2180,  
Germany national toll free 08001-536 376,  
Germany international +49 6184 90 6940, Italy +39 02 02 95059 434-254-375,  
Netherlands +31 76 571 4440, Nordic/Baltic countries +358 9 329 100,  
Russia/CIS +7 (812) 703 42 15, Spain/Portugal +34 93 223 09 18, Switzerland +41 44 454  
12 12, UK/Ireland +44 870 609 9203

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